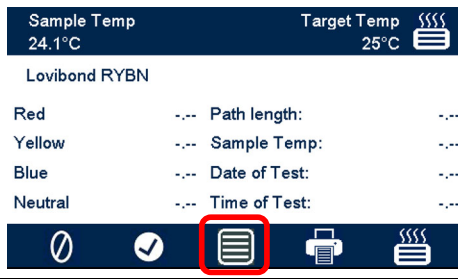
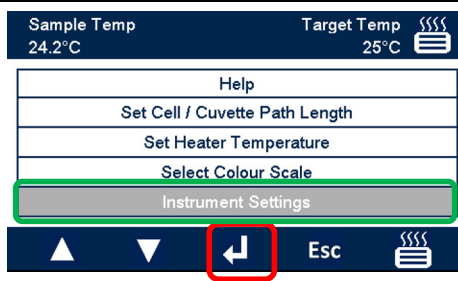

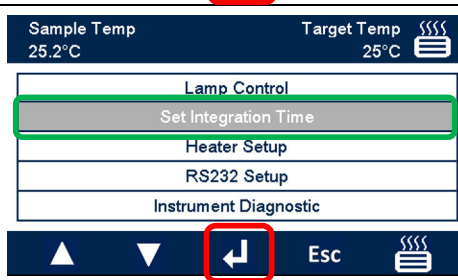
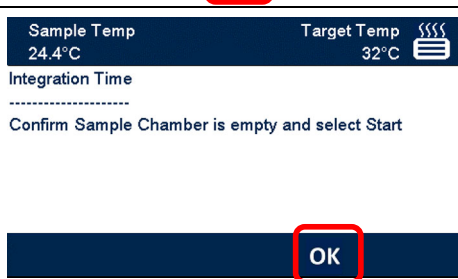
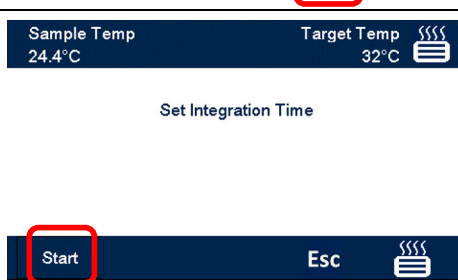


Integration Time

<p>Press the Menu Key.</p>	
<p>The Main Menu will appear. Press the Down key until "Instrument Settings" is highlighted, then press the Enter key.</p>	
<p>The Instrument Set up Menu will appear. Press the Down key until "Hardware Setup" is highlighted, then press the Enter key.</p>	
<p>The Hardware Set Up Menu will appear. Press the Down key until "Set Integration Time" is highlighted, then press the Enter key.</p>	
<p>The instrument will prompt you to check the sample chamber is empty. Press OK when confirmed</p>	
<p>Press, "Start" to begin the Integration Time process.</p>	

Time process. A progress bar will appear to show the status of the process



When the process is complete, press "OK" to return to the Hardware Set Up Menu.

